



SAFETY SHOWER AND EYE WASH STATION

Eyewash shower station is the emergency rescue facilities used in toxic and harmful environment; These devices can be emergency wash or shower when harmful chemicals splash to staff's body, face, eyes. The mainly function is to avoid further damage caused by chemical substances, but these devices can't replace medical.

CHECK ZH'S EYEWASH SHOWER STATION:

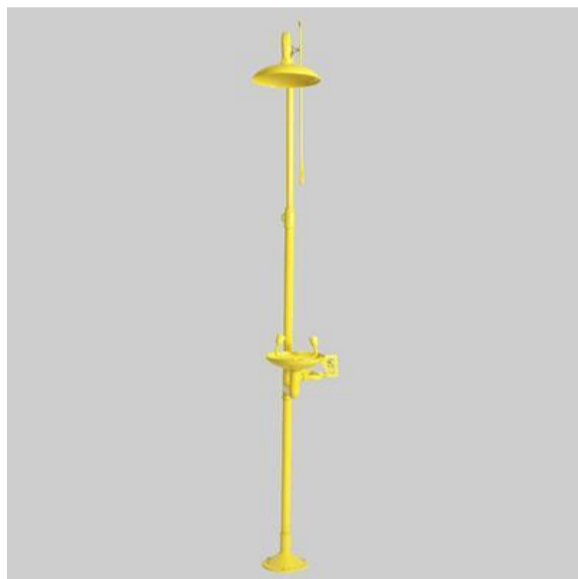

[Emergency Shower Eyewash Station](#)

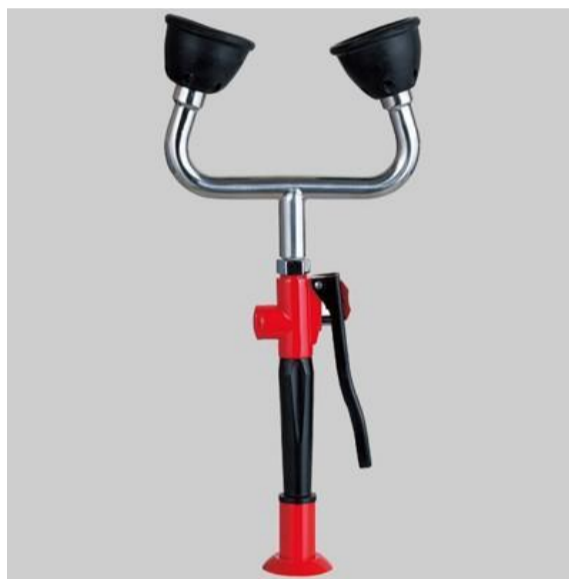

[Deck Mounted Eyewash](#)



[Portable Eyewash Station](#)

[Eyewash shower](#) is a necessary equipment for safety and labor protection, as well as an emergency and protective facility for occasions where toxic and corrosive substances such as acids, alkalis, and organic substances are exposed. There is a dust cover on each spray head of the eye washer. The valve of the eyewash is easy to operate and has anti-corrosion function. The spray head of the eye washer has a filter to prevent debris in the water from causing damage to the human eye. Meanwhile, The main valve of the emergency sprinkler device should be in a normally open state, and there should be no obstacles below the sprinkler head. It is not possible to replace the emergency sprinkler device with a normal shower device.

EYEWASH SHOWER STATION STRUCTURE

- Pedestal: used to fix the eye washer;
- Wash Area: Prevent the water from washing eyes and spraying from flowing out.
- Eyewash Shower Head: The spray system is used to spray and wash the whole body, and the eyewash system is used to wash the eyes, face, neck, or arms.
- Dust cover: A dustproof device used to protect the eye wash nozzle.
- Valve: A valve device used to open and close water flow.
- Filter: used to filter out debris entering the eye washer.

GENERAL PRINCIPLES FOR SETTING EYEWASH SHOWER STATION

- The eyewash is installed in hazardous chemical areas, and the time to reach the eyewash does not exceed 10 seconds.
- Rescue range: within 15 meters
- There are no electrical switches around the eyewash to prevent accidents.
- There should have a sewer or wastewater treatment tank.



COMBINATION EMERGENCY SHOWER AND EYEWASH STATION

We are your one-stop shop for safety [shower and eyewash station](#). ZH Lab power to honesty, high quality and relatively low cost as the base, participate in market competition, and won long-term cooperation from users.

Call now! +86-13302388256



EMERGENCY EYEWASH SHOWER STATION FEATURES

- Emergency Eyewash Shower can spray water in 1s to wash both eye and body;
- Double corrosion protection: the body is made of 304 stainless steel material coated by ABS powder with yellow finish for better chemical resistant, water resistant.
- On stock: we can ship most of in-stock items within 5 business days. Large quantity orders may require longer lead times.
- Follow Ansi z358.1



- Pedestal: aluminum alloy + yellow ABS material, high strength, and pressure resistance. Diameter 220mm, Fix with M8 expansion screw.
- Connection of inlet and outlet: DN32 internal thread, middle inlet, and lower drain. It can be customized according to requirements.
- Performance: It has good resistance to corrosive solutions such as acid, alkali, salt, oil, etc.
- Packing: 74*28*23cm;